


How to use SensiMedia for Sterility Test

MicroBio Corporation

1. General Description

This SensiMedia is designed to detect any type of microorganism as broad as possible.

2. Detection Procedure

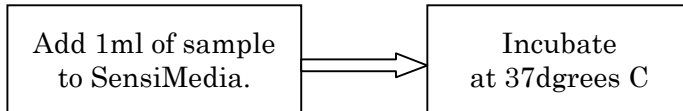

Add sample

Open a cap of SensiMedia and add 1ml of sample. Close a cap and place the SensiMedia in an incubator.

Incubate Temperature

Incubate the sample added SensiMedia at 37 degrees C.

3. Detection Criteria

The performance characteristic of liquid medium is shown in graph below. Refer to it and set up a protocol for the test. For instance, if the color of sensor remains in dark blue after 24 hours, ordinary microorganisms are not in presence (excluding mould, bacillus and the like).

