


How to use Lactose Broth SensiMedia for Coliform Detection

MicroBio Corporation

1. General Description

This SensiMedia uses Lactose Broth as liquid media.

2. Detection Procedure

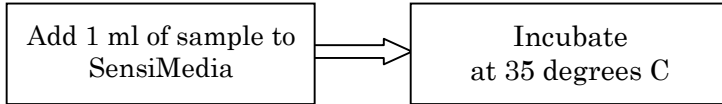

Add sample

Open a cap of SensiMedia and add 1ml of sample. Close a cap and place it in an incubator.

Incubate Temperature

Incubate the sample added SensiMedia at 35 degrees C.

3. Detection Criteria

The performance characteristic of liquid medium is shown in graph below. Refer to it and set up a protocol for the test. When the color of CO₂ sensor turns into yellowish transparent, Coliforms are detected.

