


How to use BGLB SensiMedia for Coliform Detection

MicroBio Corporation

1. General Description

This SensiMedia uses BGLB (Brilliant Green Lactose Bile) as the culture media.

2. Detection Procedure

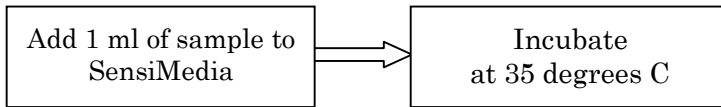

Add sample

Open a cap of SensiMedia and add 1ml of sample. Close a cap and place the SensiMedia in an incubator.

Incubate Temperature

Incubate the sample added SensiMedia at 35 degrees C.

3. Detection Criteria

The performance characteristic of BGLB is shown in the graph below. Refer to it and set up a protocol for the test. When the color of CO₂ sensor turns into yellowish transparent, Coliforms are detected.

