


How to use SensiMedia for Sterility Test

MicroBio Corporation

1. General Description

This SensiMedia is designed to detect any type of microorganism as broad as possible.

2. Detection Procedure

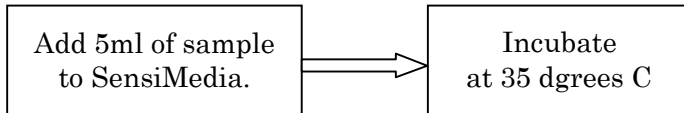

Add sample

Open a cap of SensiMedia and add 5ml of sample. Close a cap and place the SensiMedia in an incubator. If the sample is water, sample can be added to the tube as much as it is able to take.

Incubation Temperature

Incubate the sample added SensiMedia at 35 degrees C.

3. Detection Criteria

The performance characteristic of liquid medium is shown in graph below. Refer to it and set up a protocol for the test. For instance, if the color of sensor remains in dark blue after 48 hours, ordinary microorganisms are not in presence (excluding mould, bacillus and the like).

